

VIRGINIA DEPARTMENT OF FORENSIC SCIENCE

ABI 3500xl Monthly QC Log

CE _____

Task	Frequency
Use a lab wipe to clean the anode buffer container valve pin assembly on the polymer delivery pump	Monthly or as needed
Run the Wash Pump and Channels Wizard (conditioning reagent)	Monthly or as needed
Flush the pump trap	Monthly or as needed
Empty the condensation container (right of autosampler tray) and the water trap waste container (right of the pump block)	Monthly or as needed
Replace cathode buffer septa	Monthly or as needed
Clean the autosampler and drip tray	Monthly or as needed
Remove dried polymer from the capillary tips with a lint-free wipe moistened with deionized water	Monthly or as needed
Defragment the hard drive	Monthly or as needed

DATE	INITIALS	CONDITIONING REAGENT LOT#/EXPIRATION/SOURCE
		/AB

*Date/initials above indicates that ALL QC tasks for that month were completed